

SPECIFICATIONS

Range:

-58 to 302°F (-50 to 150°C)

Resolution

0.1° from -20.0 to 100°C

Accuracy:

±1°C from -20 to 200°C

OPERATION

1. Slide the power "ON/OFF" switch to the "ON" position.
2. Slide the "C/F" switch to select the desired reading.
3. Press "HOLD" button once to freeze the display.
4. Press the "HOLD" button again to resume measurement.

5. Slide the power "ON/OFF" switch to the "OFF" position to conserve battery life.
6. Keep plastic housing and plastic handle away from heat source. Do not place inside oven or expose to open flame.
7. Do not immerse probe handle or cable in liquid.

BATTERY REPLACEMENT

1. Use a coin to open the battery cover on back of unit.
2. Replace with catalog number 1039.

TRACEABLE[®]
MINI-
THERMOMETER
INSTRUCTIONS