



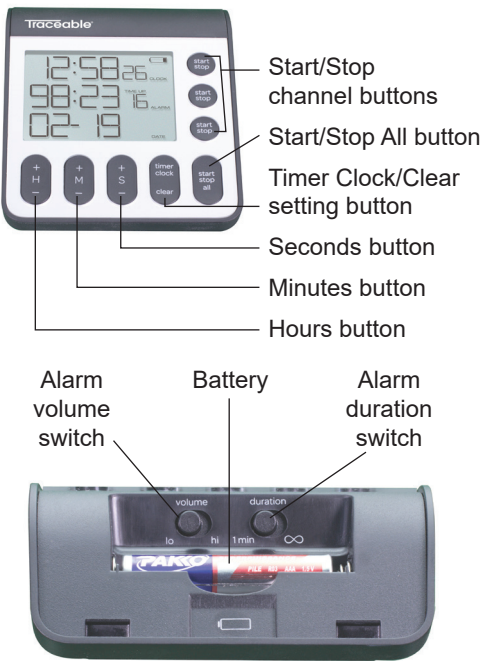


Traceable® 3-Channel Timer

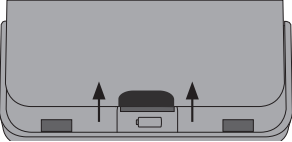


INITIAL SETUP

- 1 Remove battery cover and insert batteries.

- 2 Use the volume slider to adjust buzzer volume.

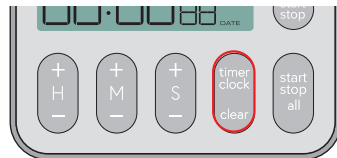- 3 Use the duration slider to adjust alarm duration.

- 4 Replace battery cover.

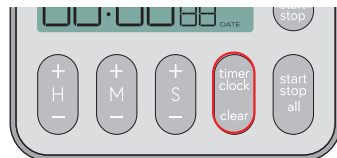
- 5 Press TIMER/CLOCK until time appears on screen.



- 6 Hold TIMER/CLOCK until there is an audible beep.



- 7 Use H/M/S buttons to set time; press TIMER/CLOCK when done.

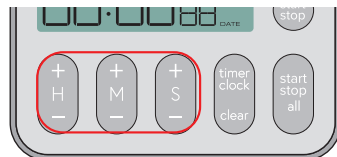


- 8 Use H/M/S buttons to configure alarm and date.



Note: To activate alarm, press START/STOP ALL button when ALARM is flashing. ALARM will appear on display if alarm is active.


SETTING A TIMER

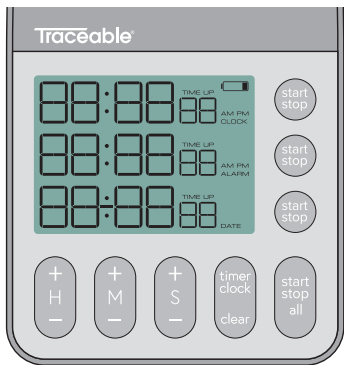

1. Press TIMER/CLOCK until timer is displayed.
2. Press and hold START/STOP on the timing channel to be configured.
3. Use H/M/S buttons to configure timer.
4. Press START/STOP on timing channel to set timer.

Note: To recall memory, simply press START/STOP button after the timer has alarmed.

STARTING A TIMER

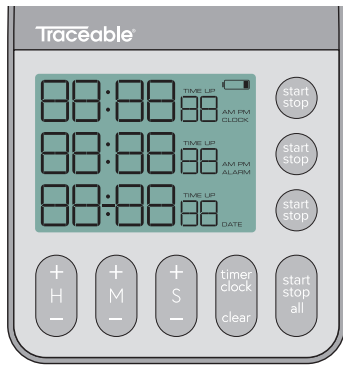

To start/stop an individual channel, press the START/STOP button for that channel.

To start/stop all channels, press the START/STOP ALL button

Note: If a channel has not been configured with a set time, pressing START/STOP on that channel or START/STOP ALL will start a count up and TIME UP will flash on that channel.

WARRANTY, SERVICE, OR RECALIBRATION

Traceable®
an antylla scientific company

12554 Old Galveston Rd. Suite B230
Webster, Texas 77598 USA
Ph. 281-482-1714 • Fax 281-482-9448
support@traceable.com
traceable.com

Traceable® Products are ISO 9001:2018 Quality-Certified by DNV and ISO/IEC 17025:2017 accredited as a Calibration Laboratory by A2LA.

Item No. 56000-13

Model No. 5665

©2022 1065T8_M_92-5665-00 Rev. 0 09022022

Traceable® 3-Channel Timer Quick-Start Guide

