sartorius

**Operating Instructions | Betriebsanleitung |  
Mode d'emploi | Instrucciones de manejo | 操作说明**

## Sartorius YRB11Z

**External Rechargeable Battery Pack | Externer Akku |  
Accumulateurs externes rechargeables | Equipo externo  
de baterías recargables | 外部可充电电池组**

98647-004-83

---

## **English      Page 3**

In cases involving questions of interpretation, the German-language version shall prevail.

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---

# Contents

<b>Contents</b> .....	3
<b>Safety Precautions</b> .....	4
Safety .....	4
<b>Operation Design</b> .....	6
Power Meter .....	6
Power Meter Explanation .....	6
<b>Installation and Operation</b> .....	7
Charging the Battery Pack .....	7
<b>Disposal</b> .....	9
<b>Specifications</b> .....	10
<b>Accessories</b> .....	10
<b>Information on Hazardous Substances</b> .....	11
<b>Declaration of Conformity</b> .....	52

---

# Safety Precautions

Please read these installation instructions carefully before putting the equipment into operation for the first time. Observe the safety instructions.

Add: The battery pack may only be used with Sartorius devices that explicitly name this battery pack YRB11Z as an accessory in their manual.

## Safety

The device complies with EU directives and standards for electrical safety and electromagnetic compatibility. Improper use or handling can, however, result in damage and/or injury. Any improper use or operation of the device, i.e., that is not consistent with the instructions, will result in forfeiture of all claims under the manufacturer's warranty.

Prevent the risk of electric short circuit.

If the battery pack becomes damaged, do not touch the battery leakage directly:

Electrolyte leakage can cause the skin discomfort. In case the electrolyte comes into contact with the eyes, rinse with water as soon as possible. Do not rub eyes. Seek medical attention as soon as possible.

### **Danger of Explosion!**

Do not use this equipment in hazardous areas in which explosive materials are present.

Make sure that the voltage rating printed on the AC adapter is identical to your local mains voltage.

---

When used in environment with electrostatic charges the connected device should be grounded.

---

Do not operate the device if the housing or AC adapter is damaged. Immediately disconnect the damaged device from the power by pulling the plug.

---

Do not expose the equipment to extreme temperatures, aggressive chemical vapors, moisture, shocks, vibration, or strong electromagnetic fields. Observe the conditions of operation described in the specifications!

---

The operator shall be solely responsible for any modifications to the equipment and for connecting any cables or equipment not supplied by Sartorius. Information on operational quality is available upon request from Sartorius. You should only use peripherals supplied by Sartorius.

---

Do not compromise the IP protection of the device.  
Do not allow liquids to enter the device.

---

Before cleaning the AC adapter or the device:  
Unplug from the mains power.

---

Lay the cables where they pose no risk of causing someone to trip.


---

Please recharge your battery pack every six months even when you do not use it for a prolonged period in order to ensure proper product performance.

---

# Operation Design

## Power Meter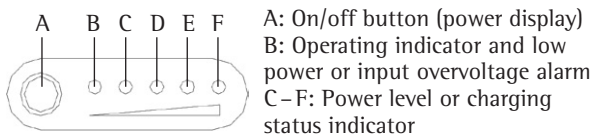

### Power Meter Explanation:

B flashing	Less than 5% power level
B flashing quickly	Input overvoltage alarm and protection
C on	About 25% power level
C+D on	About 50% power level
C+D+E on	About 75% power level
C+D+E+F on	Over 95% power level

Press the on/off button briefly: the number of LED lights will show you the power level.


---

# Installation and Operation


This battery pack is designed for use with Sartorius AC adapter GPP8 and for connection to devices that use this AC adapter. Use the AC adapter GPP8 that was supplied with your device to charge the battery pack.

For details on the AC adapter, see chapter Accessories.

## Charging the battery pack while supply power to your device (DC IN and DC OUT)

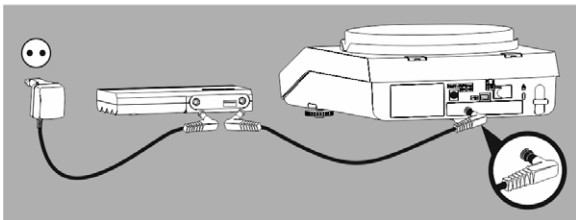
1. Connect the AC adapter to the DC IN jack of the battery pack.  
Connect the Sartorius AC adapter to the AC power source.
2. Connect one end of the DC output cable to the DC OUT jack of battery pack. Connect the other end of the DC output cable to the DC IN jack of your device.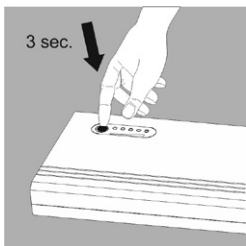


3. Press the on/off button for about 3 sec.: the battery pack will turn on.

4. The LED "B" will light up steady in red while supplying power to your device. Meanwhile, the C, D, E and F LED lights will light up in sequence while the battery pack is charging.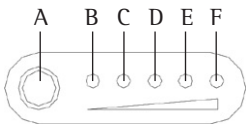

5. The C, E, E and F LED lights will light up steady when the battery pack is fully charged.

6. If AC power is not available, it will automatically switch from the AC supply source to the battery pack to supply power to your device.

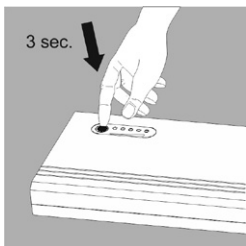7. Press the on/off button for about 3 sec.: the battery pack will turn off, and the device will be disconnected from power.

---

# Disposal

The packaging consists of environmentally friendly materials that can serve as secondary raw materials. If the packaging is no longer required, it can be disposed of in Germany at no charge via the dual system of the company VfW (contract number D-59101-2009-1129). Otherwise, dispose of the material according to the applicable local regulations for waste disposal.

The device, including accessories and batteries, should not be disposed of as household waste. It should instead be recycled as for electrical and electronic devices.

For more information regarding disposal and recycling, please contact our local service representatives.

In addition, the partners listed on the following website are available to assist you within the EU:

- 1) Load <http://www.sartorius.com>.
- 2) In the header, type "Service".
- 3) Then select "Disposal Information".
- 4) The addresses of the local Sartorius disposal contact persons can be found in the PDF files listed on this page.

Devices contaminated with hazardous materials (NBC contamination) will not be accepted for repair or disposal. Detailed information with service addresses for returning your device for repair or disposal can be found on our website ([www.sartorius.com](http://www.sartorius.com)) or requested from a Sartorius Service Center.

---

## Specifications

Model	YRB11Z
Capacity	49 Wh (4.4 Ah/11.1 V)
Input	DC 15 V, 500 mA
Output	DC 15 V/1 A, 5 V/1 A
Standby Current	< 25 $\mu$ A
Operating Temperature	0 to +40°C
Storage Temperature	-10 to +45°C
Operation Humidity	$\leq$ 95%
Storage Humidity	45% – 75%
Size	135 mm $\times$ 77 mm $\times$ 25 mm
Net Weight, approx.	380 g
Packaging Weight	500 g
Charging Time with GPP8	8 hours (20°C)

## Accessories

AC adapter for the battery charger GPP8:

- AC adapter versions Europe, USA, Australia, India, South Africa 6971790
- AC adapter versions Korea, Brazil, Argentina 6971791

---

# Information on Hazardous Substances

<b>UN number</b>	<b>Road Transport in Europe ADR</b>	<b>International Marine Transport incl. RoRo Transport IMDG Code</b>	<b>Air Transport International IATA-DGR</b>
<b>UN 3480</b>	SV 188	SV 188	PI 965
Lithium ion battery	SV 230	SV 230	
Max. 20 Wh/cell Max. 100 Wh/batt.	“Simplifying” rules	“Simplifying” rules	“Simplifying” rules

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*This declaration certifies conformity with the above mentioned EU Directives, but does not guarantee product attributes. Unauthorised product modifications make this declaration invalid. The safety information in the associated product documentation must be observed.*

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Sartorius Lab Instruments GmbH & Co. KG  
Goettingen, 20/04/2016

Dr Reinhard Baumfalk  
Vice President R&D

Dr Dieter Klausgrete  
Head of International Certification Management

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Sartorius Lab Instruments GmbH & Co. KG  
Goettingen, 20/04/2016

Dr. Reinhard Baumfalk  
Vice President R&D

Dr. Dieter Klausgrete  
Head of International Certification Management

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Phone +49.551.308.0  
Fax +49.551.308.3289  
www.sartorius.com

Sartorius Scientific Instruments (Beijing) Co.,  
Ltd.No. 33 Yu An Road,  
Shunyi District, Zone B,  
Tianzhu Airport Industry Park,  
101300 Beijing, P.R. China  
Phone +86.10.8042.6300  
www.sartorius.com.cn

Copyright by Sartorius, Goettingen, Germany.  
All rights reserved. No part of this publication may  
be reprinted or translated in any form or by any means  
without the prior written permission of Sartorius.  
The status of the information, specifications and  
illustrations in this manual is indicated by the date given  
below. Sartorius reserves the right to make changes to  
the technology, features, specifications and design of the  
equipment without notice.

Status:

April 2016, Sartorius Lab Instruments GmbH & Co. KG,  
Goettingen, Germany