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## Warranty

In accordance with IKA warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

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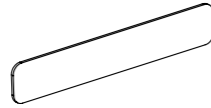
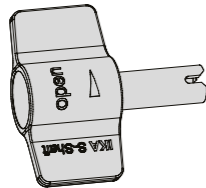
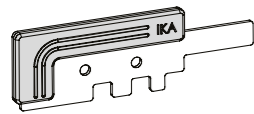
Flachschlüssel 25 F \*  
Flat key 25 F \*  
Clé plate 25 F \*

\* Nur für S 25 KV - 25 F!  
\* Only for S 25 KV - 25 F!  
\* Uniquement pour S 25 KV - 25 F!

Fig. 1

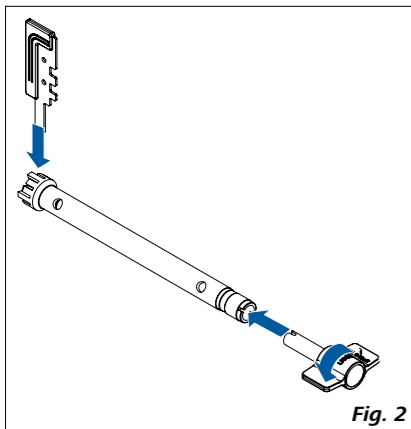
Fig. 2

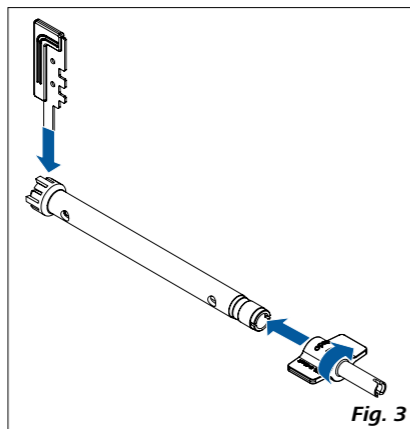
Fig. 3

# IKA

designed for scientists

### IKA-Werke GmbH & Co. KG

Janke & Kunkel-Straße 10, 79219 Staufen, Germany  
Phone: +49 7633 831-0, Fax: +49 7633 831-98  
eMail: sales@ika.de

#### USA

IKA Works, Inc.  
Phone: +1 910 452-7059  
eMail: usa@ika.net

#### MALAYSIA

IKA Works (Asia) Sdn Bhd  
Phone: +60 3 6099-5666  
eMail: sales.lab@ika.my

#### JAPAN

IKA Japan K.K.  
Phone: +81 6 6730 6781  
eMail: info\_japan@ika.ne.jp

#### VIETNAM

IKA Vietnam Company Limited  
Phone: +84 28 38202142  
eMail: sales.lab-vietnam@ika.com

#### KOREA

IKA Korea Ltd.  
Phone: +82 2 2136 6800  
eMail: sales-lab@ika.kr

#### CHINA

IKA Works Guangzhou  
Phone: +86 20 8222 6771  
eMail: info@ika.cn

#### INDIA

IKA India Private Limited  
Phone: +91 80 26253 900  
eMail: info@ika.in

#### BRAZIL

IKA Brazil  
Phone: +55 19 3772 9600  
eMail: sales@ika.net.br

#### POLAND

IKA Poland Sp. z o.o.  
Phone: +48 22 201 99 79  
eMail: sales.poland@ika.com

#### UNITED KINGDOM

IKA England LTD.  
Phone: +44 1865 986 162  
eMail: sales.England@ika.com

Discover and order the fascinating products of IKA online:  
[www.ika.com](http://www.ika.com)

IKAworldwide

IKAworldwide /// #lookattheblue

@IKAworldwide

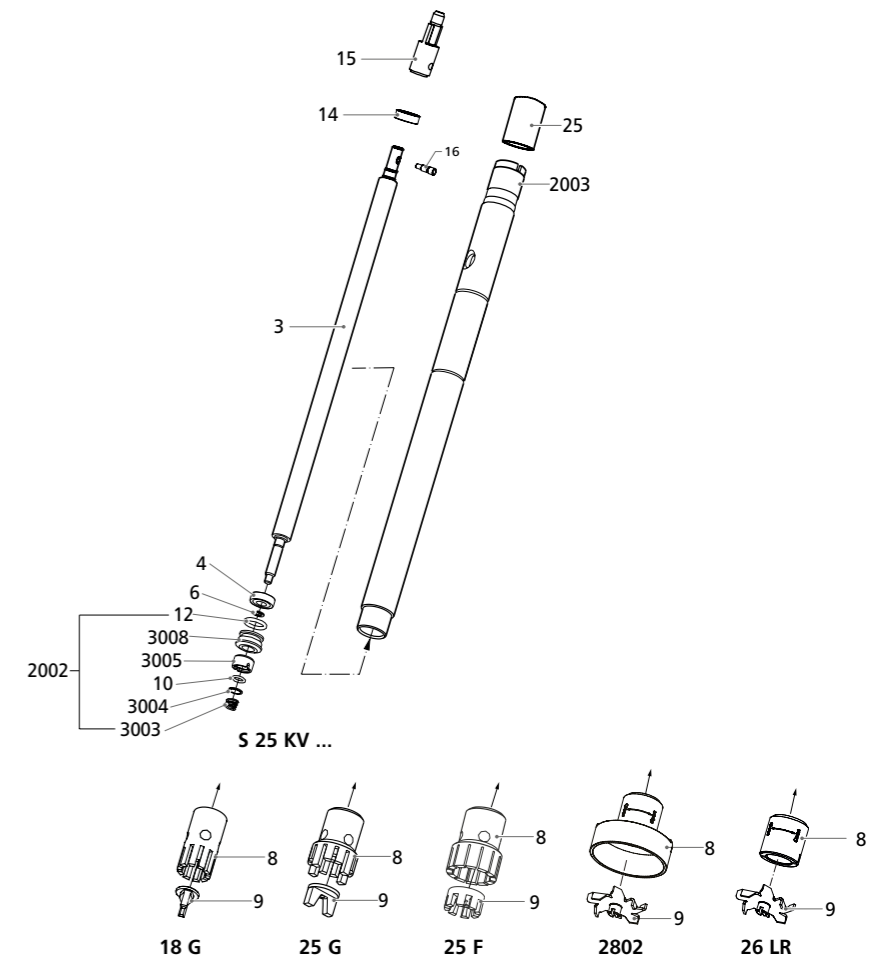
Technical specifications may be changed without prior notice.

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Source language: German

### Safety instructions

• **Read the operating instructions in full before starting up and follow the safety instructions.**

• Keep the operating instructions in a place where they can be accessed by everyone.

• Ensure that only trained staff work with the appliance.

• Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

- The dispersion tool may heats up during operation. In particular in the processing of hot medium, the tool must be cooled before it can be removed from the drive.
- When in operation the dispersion tool can become hot. If the dispersion tool is not inserted into the drive flange correctly or the medium is hot, it can become extremely hot.

• Don't touch rotating parts during operation.

• No liquid may get into the drive unit of the disperser.

• The work has to be interrupted immediately if you notice unusual noise and/ or emission of liquid from the upper rinsing drilling. The seals must be controlled, and replaced if necessary.

• While working with the disperser, the user must wear his personal protective equipment, in particular safety goggles.

• Follow the health and safety regulations accident prevention regulations.

• Note the operating instructions of the disperser unit.

• Abrasion of the dispersion equipment or the rotating accessories can get into the medium you are working on.

• **Attention!** Never run the dispersion tool dry. Without cooling by the medium the bearing will be destroyed.

• The dispersion tool is not suitable for continuous operation.

• Dispersing causes the medium to heat up.

• Likewise the optimal dispersion duration and rotating frequency must be determined by attempts.

• The function of the dispersion tools depends on the condition of the sharp edges of rotor and stator. Abrasive media can round off these edges fast, whereby the disperse effect diminishes.

#### Product information

The shaft bearing of the dispersion tool consists of a two ball bearings and a single-thrust slide ring seal (SiC).

Content of package: Dispersion tool and tool kit.

The plastic materials used are approved by FDA (Food and Drug Administration Approval).

#### Application instructions

**Application:** Wet crushing, suspension, emulsifying (Batch-operation).

#### Spare parts list

<b>Item</b>	<b>Designation</b>	<b>Item</b>	<b>Designation</b>	<b>Item</b>	<b>Designation</b>
<b>3</b>	Rotor shaft	<b>12</b>	O-Ring 9.25 x 1.78	<b>2003</b>	Shaft tube, compl.
<b>4</b>	Deep-groove ball bearing	<b>14</b>	Deep-groove ball bearing	<b>3003</b>	Pressure spring, right
<b>6</b>	Seeger circlip ring	<b>15</b>	Coupling	<b>3004</b>	Thrust collar
<b>8*</b>	Stator	<b>16</b>	Threaded pin	<b>3005</b>	Slide ring
<b>9*</b>	Rotor	<b>25</b>	Plastic cap	<b>3008</b>	Counter fit mounting, compl.
<b>10</b>	O-Ring 4.7 x 1.42	<b>2002</b>	Friction ring seal, compl.		

für **S 25 KV 2802: 8\*** – Protection ring, **9\*** – Dissolver

#### Maintenance and cleaning

Allowed procedures	Methods of sterilisation	Notice
Chemical procedures	By germ - killing dissolvents alcohol, phenol,formalin.... Disinfectant remainders are to be removed with germ-free water.	For ball bearings <b>not</b> allowed. Note the chemical resistance of ball bearings and o-rings.
Hot air	Germ - killing by hot-air with 160 to 190°C (approx. 30 min)	Life span of the bearings shortens.

For pre-cleaning the dispersion tool, operate it in in a dissolvent for dissolving the residuaries of the substance. Due to the large flow rate, rotor and stator are cleaned to a large extent. The dispersion tool must be divided and cleaned immediately after working.

For cleaning purposes disassemble the dispersion tool as described below (items see front page):

• Unscrew the rotor **(9)** from the shaft **(3)**, see **Fig. 2**:

Hold the rotor **(9)** with the flat key. Screw off the rotor **(9)** by turning the shaft **(3)** in counterclockwise using the shaft key.

• Unscrew the stator **(8)** from the shaft tube **(2003)** (left-hand thread), see **Fig. 3**:

Hold the stator with the flat key. Screw off the stator by turning the shaft tube **(2003)** in clockwise using the shaft key.

In doing so, the dispersion tool and the stator should be held up in vertical position, thus preventing the spring of the slide ring sealing from falling out.

• Tap the shaft **(3)** and the ball bearing **(4)** downward in the direction of the stator thread out of the shaft tube.

• In this position the parts of the slide ring seal can be removed from the shaft **(3)**.

• The slide rings and O-rings must be replaced in the case of wear. The sliding surfaces of the slide and counter ring must be clean.

The assembly of the dispersion tool takes place in reverse order. Make certain that the O-ring **(10)** and the thrust collar **(3004)** are properly fitted during assembly.

**Attention!** The rotor must not be tightened too strongly, because otherwise the thread will be damaged.

#### Technical Data

		<b>S 25 KV - 18 G</b>	<b>S 25 KV - 25 G</b>	<b>S 25 KV - 25 F</b>	<b>S 25 KV 2802</b>	<b>S 25 KV - 26 LR</b>
Working range	<b>l</b>	0.01–1.5	0.05–2	0.1–2	0.1–2	0.1–2
Stator/Rotor (Protection ring/Dissolver) (Ø)	<b>mm</b>	18 / 12.7	25 / 17	25 / 18	39 / 30	20 / 30
Gap between rotor and stator	<b>mm</b>	0.3	0.5	0.5	-	-
Max. allowable speed	<b>rpm</b>			25000		
Max. circumferential speed	<b>m/s</b>	16.6	22.2	23.6	39.3	39.3
Min./max. immersion depth	<b>mm</b>			40 / 225		
Material in contact with medium				FFPM / SiC, AISI 316L		
pH range				2–13		
Suitable for solvents				Yes		
Suitable for abrasive substances				No		
Max. working temperature	<b>°C</b>			220		
Sterilization methods				all meththods		
Working range vacuum	<b>mbar</b>			1		
Working range pressure	<b>bar</b>			6		
Ultimate fineness, suspensions	<b>µm</b>	10–50	15–50	5–25	-	-
Ultimate fineness, emulsions	<b>µm</b>	1–10	1–10	1–5	-	-

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