Refracto – reliable refractive index wherever you need it

Refracto 30PX and Refracto 30GS are compact, easy-to-use refractometers for measurements in laboratory and production environments. Depending on the application, these instruments can be used in two ways: Either put it on a flat surface and place a drop of sample onto the measurement cell, or immerse the sensor directly into the fluid. As soon as you press the measurement key, the result is displayed in the desired units on the backlit LC-Display.

You no longer need to rely on error-prone readings of dark/ bright transitions as with optical instruments. Refracto even measures dark samples correctly. Because the refractive index is temperature dependent, Refracto also automatically compensates the result.

The Refracto 30PX has a measuring cell made of optical glass, whereas the Refracto 30GS has a measuring cell made of sapphire. Sapphire has a higher refractive index and a better thermal conductivity than glass. For this reason the Refracto 30GS has an extended measuring range and registers the sample temperature more quickly.

PortableLab™ – Lab power in your hands! Refracto is delivered in a special case which contains everything you need for successful measurements: Pipettes for sample handling, cleaning towels, flasks for samples and cleaning solutions, operating instructions and a CD-ROM with data transmission software, a tutorial to







Good-bye Abbé: Easy-to-perform refractive index measurements



A refractive index determination is the simplest way to quickly determine the quality of a substance. Using Refracto, the determination is faster and less prone to errors than traditional methods (e.g. Abbé-refractometer). Refracto can also store, print and transfer results to a PC.







Reliable quality control



Wherever syrup concentrates, saline solutions or vinegar are produced: Refractometers are essential tools for auality control. The refractive index helps you make sure everything has gone right in the production process. With Refracto, you can perform your quality checks directly in the production facility.



Grapes and fruit: Direct display

With Refracto, you can not only display sugar content in Brix%, but also choose direct conversion into Oechsle, T.A. 1990, KMW, Baume HFCS42 and HFCS55, which are shown directly after the measurement



Quick concentration determination

The measurement of the refractive index is the quickest and easiest method to check antifreeze concentration. Refracto will tell you within seconds whether the antifreeze contains the right amount of ethylene glycol, propylene glycol, ethanol or sodium chloride. It even allows you to directly read the freezing point of these fluids in °C or °F.



Soft drinks:

Simple sugar content determination



The precise measurement of the temperature during measurement and the automatic temperature compensation using built-in ICUMSA tables ensure highly reliable Brix% results.





With Refracto, any concentration measurement can be performed without manual calculation. Just define the corresponding formula and the instrument gives you the

Refracto: Desktop features in a handheld instrument

METTLER TOLEDO

Plain language interface

Measurement cell with temperature sensor

Refracto determines refractive index using the total reflection

method. The glass prism of the measurement cell of the Re-

fracto 30PX (right) is held by a stainless steel ring, whereas

ring of hard gold plated brass. Sapphire has a higher refrac-

tive index and a better thermal conductivity than glass. For

this reason the Refracto 30GS has an extended measuring

range and registers the sample temperature more quickly.

the prism of the Refracto 30GS (left) is made of sapphire, the

The backlit LC-Display of Refracto shows your results

and settings in plain language. The simple and intuitive

menu structure allows you to change the settings within

Save and transfer your data

Refracto stores up to 1100 results

including sample identification, mea-

surement unit, temperature compensa-

tion coefficient, date and time. You are

free to transfer the data (together with

whenever you want

an instrument iden-

tification) to a PC or printer any time using the infrared interface. The PC

software to do

this comes with

Refracto.

seconds. You hardly need an instruction manual.

Ease-of-use

The clearly labelled keyboard makes operation easy and efficient

Temperature compensation

The refractive index of a sample depends on temperature. During measurement, Refracto determines the temperature and then corrects the refractive index to a standard temperature of 20 °C or any other temperature the user defines. To make quick measurements of different types of samples, you can easily switch between up to 10 user defined correction coefficients

It is your choice: hand-held or bench-top

Refracto is a hand-held and a benchtop refractometer in one. You either place the instrument on a flat surface and add a drop of sample onto the measurement cell or immerse the cell directly into the sample.

PortableLab™

Specifications Refracto 30PX / Refracto 30GS

Measurement method: Determination of the angle of total reflection of the D-line of sodium (589.3 nm) • Refractive index: Measurement range: 1.32 - 1.50 (PX), 1.32 – 1.65 (GS), Resolution: 0.0001, Accuracy: ±0.0005 • Brix%: Measurement range 0 – 85 Brix%, Resolution: 0.1 Brix%, Accuracy ± 0.2 Brix% • Temperature: Measurement range 0 – 85 Brix% Resolution: 0.1 Brix%, Accuracy: ±0.0005 • Brix% • Temperature: Measurement range 0 – 85 Brix% Resolution: 0.1 Brix%, Accuracy: ±0.2 Brix% • Temperature: Measurement range 0 – 85 Brix% Resolution: 0.1 Brix%, Accuracy: ±0.0005 • Brix% • Temperature: Measurement range 0 – 85 Brix% • Temperature: ment range: 10 - 40 °C, Resolution: 0.1 °C, Display: °C or °F, Ambient temperature: 5 - 35 °C • Units of measurement: nD, nD temperature compensated, Brix%, HFCS42, HFCS55, T.A. 1990, "KMW (Babo), "Baume, "Oechsle (D, CH), w/w%, v/v%, spec. gravity and freezing point ("C or "F) for ethanol and NaCl, w/w%, v/v% and freezing point (°C or °F) for ethylene glycol and propylene glycol, w/w% and v/v% for isopropanol, user defined unit • Temperature compensation: With user-defined temperature compensation coefficient (nD temperature compensated, user defined), or automatically (all other units). Up to 10 temperature correction coefficients can be stored in the instrument • Calibration: With pure water • Data memory: For up to 1100 results (result with unit, sample identification, temperature correction coefficient, date and time) • Display: Backlift LC-Display • Interface: Infrared for data transfer to PC and printer (IrDA or RS232C) • Weight: Approx. 250 a • Batteries 2 x LR3, 1.5 V, type AAA approx. 60 hours battery life (one measurement per minute) • Materials: Housing: PBT. Measurement cell (PX): Glass, stainless steel. Materials with sample contact (PX): Glass, stainless steel, PBT, Measurement cell (GS): Sapphire, hard gold plated brass, Materials with sample contact (GS): Sapphire, gold, PBT,

PortableLab"



Refracto 30GS

www.mt.com/refracto

portable hiah-end refractometer

Refracto 30PX www.mt.com/refracto

portable refractometer

Densito 30PX

www.mt.com/densito

density meter

Seven2Go

www.mt.com/Seven2Go

Portable instruments

for pH, lon, Cond, and DO measurements



Quality certificate. Development, production and testing according to ISO 9001. Environmental management system according to ISO 14001.



Worldwide service. Our extensive service network is among the best in the world and ensures maximum availability and service life of your



European conformity. The CE conformity mark provides you with the assurance that our products comply with the most recent EU directives.



On the Internet. You will quickly find lots of essential information about our products, our services, and our company at

www.mt.com

METTLER TOLEDO

gone global...

the contact addresses of METTLER TOLEDO representatives globaly can be found under the Internet address www.mt.com/contacts

otherwise:

Mettler-Toledo GmbH

PO Box VI-400 CH-8606 Greifensee Phone +41-44-944 22 11, Fax +41-44-944 31 70

Mettler-Toledo GmbH, Analytical

CH-8603 Schwerzenbach, Switzerland Phone +41 44 806 77 11 Fax +41 44 806 73 50 Internet: www.mt.com

Subject to technical changes © 12/2014 Mettler-Toledo GmbH, 30243620

Refracto 30PX / 30GS Refractometer Refracto METTLER TOLEDO



