


## Thermo Scientific Nalgene Forceps

Cat. No. 6320-0010 (blue polypropylene)

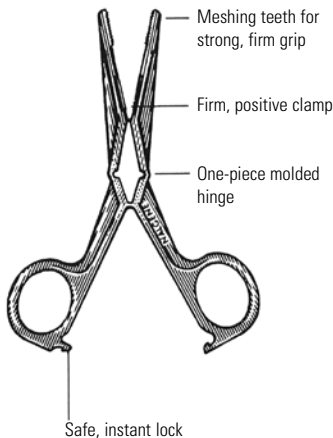

Disposable or reusable.  
Autoclavable up to 121°C/15 psig;  
can also be sterilized by ethylene  
oxide gas.

**IMPORTANT:** Sterilize and store  
forceps **IN THE OPEN POSITION** to  
preserve the natural tension of the  
one-piece hinge.

Nalgene® Forceps are useful for  
handling Nalgene Cryogenic Vials,  
Cat. No. 5000. They can also be  
used as clamps for  
chromatography columns.