





»I keep cool in tough situations, so do my samples.«

The new Centrifuge 5425 R – Your Lifelong Companion



Your Lifelong Companion

Refrigerated Centrifuge 5425 R is the new laboratory standard for temperature sensitive samples up to 5 mL. It features state of the art cooling technology to ensure 4 °C at full speed for maximum sample protection and best separation results. Our solution for your everyday challenges.

Benefits

- > More applications: Rotor portfolio extended from 4 to 6 including 5 mL tube rotor and swing-bucket rotor for PCR strips
- > More flexibility: Rotors interchangeable between air-cooled Centrifuge 5425 and new refrigerated Centrifuge 5425 R
- > More throughput: Maximum sample volume per tube increased from 2 mL to 5 mL



keypad control

version

- rotary knob version
- Quick setting of parameters or easy cleaning of keys

> More versatility:

Available with rotary knob

and keypad control -

New Features



Enhanced convenience

- > One push short-spin key no need to hold key all the while
- > End of run timer (time after rotor stopped)



Improved ergonomics

- > Eppendorf QuickLock[®] rotors for ergonomic lid locking
- > Closes with only a ¼ turn and takes repetitive stress off your wrist



Increased versatility

- > 10-place 5.0 mL rotor for screw cap and snap lid tubes
- > Swing-bucket rotor for divisible plates to speed up your PCR/qPCR setup



5 mL sample volume

Special feature: Accommodates adapters to hold 1.5/2 mL, HLPC, and Cryo tubes Aerosol-tight: For safe centrifugation of hazardous samples Max. capacity: 10 × 5 mL tubes (screw cap & snap lid) Max. speed: 21,300 × g (15,060 rpm)

Rotor FA-10×5



Rotor S-96×0.2-PCR

Special feature: Swing-bucket rotor: Ideal for small volume PCR/qPCR experiments as samples are collected in the center of tubes **Max. capacity:** 96 x PCR tubes with 0.2 mL or $4 \times \frac{1}{4}$ divisible Eppendorf twin.tec[®] PCR plate 96 (unskirted) **Max. speed:** 3,217 × *g* (6,000 rpm)

Spin column based DNA & RNA purification with tube lids open

Unique feature: Designed with an extra high rim to support open tube lids during centrifugation Aerosol-tight: For safe centrifugation of hazardous samples Max. capacity: $18 \times 1.5/2.0$ mL tubes Max. speed: $18,565 \times g$ (15,060 rpm)

> Micro-volume protocols, such as PCR/ qPCR reaction set-up in 0.2 mL tubes and strips

Special feature: Windshield design for reduced turbulences during the run for whisper quiet operation. **Max. capacity:** $32 \times PCR$ tubes or $4 \times PCR$ strips with 0.2 mL **Max. speed:** $18,134 \times q$ (15,060 rpm)

Pre/Post-PCR and qPCR processing with up to 96 samples per run

Centrifuge Profile



Enhanced convenience Soft-touch one-finger lid closure for ergonomic

Reduced noise pollution

OptiBowl[®] design for quiet operation, even without rotor lid

Longer lasting

Stainless steel rotor chamber with built-in condensation drain to eliminate water accumulation

Superior temperature control

Aluminum rotors, with high heat conductivity for fast pre-cooling

Improved comfort

New easy to read super bright display

eppendorf

<u> </u>		
Ordering	intorm	nation
oracing	morn	ation

Description	Order no.
Centrifuge 5425 R, keypad control, with rotor FA-24×2, 230 V/50-60 Hz	5406 000 119
Centrifuge 5425 R, rotary knob control, with rotor FA-24×2, 230 V/50-60 Hz	5406 000 518
Rotor	
Rotor FA-24x2	5495 500 006
Rotor FA-24x2-PTFE	5495 503 005
Rotor FA-10x5	5495 505 008
Rotor FA-18×2-Kit	5495 508 007
Rotor F-32×0.2-PCR	5495 510 001
Rotor S-96×0.2-PCR	5495 512 004
Adapter	
Adapter, for 1 tube 0.5 mL or microtainer, max. Ø 8 mm, for all 1.5/2.0 mL rotors, set of 6 pcs.	5425 716 001
Adapter, for 1 PCR tube 0.2 mL, max. Ø 6 mm, for all 1.5/2.0 mL rotors, set of 6 pcs.	5425 715 005
Adapter, for 1 vessel 1.5-2.0 mL, for all 5.0 mL rotors, set of 4 pcs.	5820 768 002
Adapter, for 1 cryo tube, for all 5.0 mL rotors, set of 4 pcs.	5820 769 009
Adapter, for 1 HPLC tube, for all 5.0 mL rotors, set of 4 pcs.	5820 770 007
Adapter, for 1 tube 0.4 mL, max. Ø 6 mm, for all 1.5/2.0 mL rotors, set of 6 pcs.	5425 717 008

Premium Services for Your Centrifuges



Centrifuge Performance Plans

An annual inspection and maintenance of centrifuges and rotors by an Eppendorf certified service technician is highly recommended to ensure a safe working environment.

Certification Services

Installation Qualification (IQ) and Operational Qualification (OQ) services make sure that temperature accuracy, rotational speed verification and operating time meet manufacturer's specifications.

Further information and local offers in your country.



www.eppendorf.com/epServices



Your local distributor: www.eppendorf.com/contact Eppendorf AG · Barkhausenweg 1 · 22339 Hamburg · Germany eppendorf@eppendorf.com · www.eppendorf.com

www.eppendorf.com/cool-in-tough-situations

Eppendorf[®], the Eppendorf Brand Design, the epServices logo, Eppendorf QuickLock[®], OptiBowl[®] and Eppendorf twin.tec[®] are registered trademarks of Eppendorf AG, Germany. U. S. Design Patents are listed on www.eppendorf.com/ip. All rights reserved, including graphics and images. • Copyright © 2020 by Eppendorf AG. Order No.: AA01 014 320/EN1/GLP/PDF/SSO