

INTEGRA



PIPETBOY pro Enlightened pipetting

PIPETBOY pro


Liquid handling lies at the core of every life science experiment and a key to successful research is the ability to handle liquids with precision and efficiency.

The PIPETBOY pro pipetting aid unites all essential characteristics for productive work with serological pipettes: accuracy, speed and ergonomic design.

Accurate and fast filling of all pipette types

The accuracy of pipetting depends on precise graduations on the serological pipette and on the ability to accurately match the liquid's meniscus with the graduation mark. The PIPETBOY pro enables the user to very accurately displace liquid in the pipette by applying variable finger pressure on the pipettor triggers.

Fully press the aspiration trigger to quickly draw liquid into the pipette until almost full, then gradually release the trigger to reduce speed and accurately match the meniscus to the graduation mark. Similarly, when dispensing, slightly press the trigger for drop-by-drop dispensing of liquid (gravity dispensing), or fully trigger the powerful pump motor for fast emptying of the pipette.

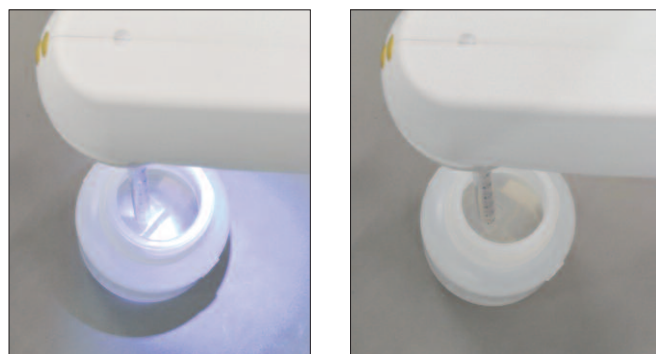

The PIPETBOY pro with the LED switched on (left) makes the bottom of the bottle clearly visible compared to the bottle that is not illuminated (right).



Illumination of the work area

Adequate lighting is a prerequisite for error-free and comfortable pipetting. The LED integrated in the PIPETBOY pro provides an optimal illumination of the area around the pipette, ensuring accurate pipetting and preventing eyestrain. The LED lights up only when pressing the pipetting triggers and automatically switches off after a few seconds. You can also turn off the illumination function completely if necessary (Patent pending).

Quality you can rely on

PIPETBOY pipette controllers are designed and manufactured to the highest quality standards and therefore offered with an extended 3 year warranty upon registration.



Design matters

The PIPETBOY pro has been designed following the recommendations of hand shape ergonomics. The result is an extra-light pipettor of 190 grams that comfortably fits in your hand and makes even prolonged pipetting effortless.

To guarantee a firm grip of small and large pipettes alike, the pipette mount in the nose of the PIPETBOY pro has a special conical shape which can be inverted when a one or two ml pipette is used creating perfect stability.

The PIPETBOY pro is available in five colors so that you can assign a specific color to a particular application and so avoid mix-ups. You can also simply pick your favorite one and brighten your lab.

The PIPETBOY pro is powered by high-quality NiMH batteries (standard AAA type) that you can charge approximately one thousand times. When fully charged, the PIPETBOY pro can accomplish 4000 pipetting cycles of 25 ml each before it needs recharging. Once the charge becomes nearly depleted, a red indicator light alerts you when batteries need to be charged.




Smart accessories

Place your PIPETBOY pro in the **charging station** when not in use and the batteries will be automatically charged, so that you always have a pipettor that is ready to use.

The **stand** for the PIPETBOY and VACUBOY is the ideal accessory for working in a cell culture hood. It keeps the workplace organized and the PIPETBOY pro, with or without inserted pipette, is always at hand.

Stable positioning

The smart design of PIPETBOY pro allows you to place the unit stably on its back side, even with a serological pipette inserted.

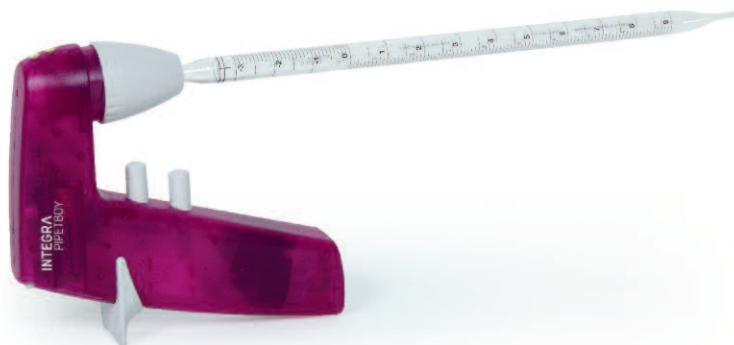

Technical data

Pipetting specifications

Maximum pipetting speed	12 ml/s
Presetting of speed (via thumb wheel)	2 – 12 ml/s
Pipette compatibility	glass and plastic pipettes with volumes from 1 to 100 µl
Filter protection	0.45 µm (included) or 0.2 µm (order separately)
Weight	190 g
Gravity dispensing	slight pressure on dispense button

Batteries charging specifications


Autonomy (of fully charged batteries)	4000 aspiration and dispensing cycles of 25 µl.
Charging of NiMH batteries	Up to 1000 times per battery. User can replace batteries.
Charge indicator light	Red: charging needed Green: connected to current
Rechargeable battery	Widely available AAA battery (1.2V, 1000 mAh)

Ordering information

Instrument		Theme	Part No.
PIPETBOY pro	Pipettor complete with spare sterile filter, two batteries, mains adapter and wall mount.	white	156 400
		pink	156 401
		green	156 402
		orange	156 403
		blue	156 404

Accessories		Part No.
Charging station	for parking and charging the PIPETBOY pro	156 450
Stand	for PIPETBOY and VACUBOY	155 065
Mains adapter (100-240 VAC, 50/60 Hz)	US version	156 630
	EU version	156 631
	UK version	156 632
	AU version	156 633
	JP version	156 634

Consumables		Part No.
Filter 0.45 µm	sterile	153 015
	non-sterile	153 016
Filter 0.2 µm	sterile	156 608
	non-sterile	156 607
Pipette mount	for holding the pipette in the nose piece, silicone	151 020
Battery set	2 AAA batteries, rechargeable	156 151

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

INTEGRA Biosciences AG
7205 Zizers, Switzerland
T +41 81 286 95 30
F +41 81 286 95 33
info@integra-biosciences.com

INTEGRA Biosciences Corp.
Hudson, NH 03051, USA
T +1 603 578 5800
F +1 603 577 5529
info-us@integra-biosciences.com

INTEGRA Biosciences Ltd.
Egham, Surrey TW20 9EY, UK
info-uk@integra-biosciences.com



www.integra-biosciences.com

INTEGRA