20000005114

**IKA**<sup>®</sup>

UTTD 122014

## IKA<sup>®</sup> ULTRA -TURRAX<sup>®</sup> Tube Drive

Betriebsanleitung Ursprungssprache	DE	4
Operating instructions Mode d'emploi	EN FR	11 18
Руководство по эксплуатации	RU	25
Indicaciones de seguridad Veiligheidsinstructies Norme di sicurezza Säkerhetsanvisningar Sikkerhedshenvisninger Sikkerhetsanvisninger Turvallisuusohjeet Instruções de serviço Wskaźówki bezpieczeństwa Bezpečnostní pokyny Biztonsági utasítások Varnostna opozorila Bezpečnostné pokyny Ohutusjuhised Drošības norādes Nurodymai dėl saugumo	ES NL IT SV DA NO FI PT PL CS HU SL SK ET LV LT	32 33 34 35 36 37 38 39 40 41 42 43 44 45 46 47
Указания за безопасност Instrucțiuni de securitate Υποδείξεις ασφαλείας	BG RO EL	48 49 50

IKA-Works, Germany  
Reg. No. 004343

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### 1 EC-Declaration of conformity

We declare under our sole responsibility that this product corresponds to the regulations 89/336EC, 98/37EC and 2006/95EC and conforms with the standards or standardized documents DIN EN IEC 61010-1, DIN EN IEC 61326-1, DIN EN ISO 12100-1, -2 and EN 60204-1.

### 2 Explanation of symbols

General hazard

This symbol identifies information **that is of vital importance for safeguarding your health and safety**. Disregarding this information can lead to health impairment and injuries.

This symbol identifies information **that is of importance for the technically correct functioning of the system**. Disregarding this information can result in damage to the instrument or to system components.


Risk of biohazard infection.

Risk of hand injury.

### 3 Safety instructions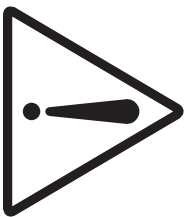

*For your protection*

#### • Read the operating instructions in full before starting up and follow the safety instructions.

- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- The instrument must be supervised at all times when in operation.

- **Caution!** For research purposes only! Not suitable for diagnostic use (in accordance with the IVD Directive).

**DANGER**

The instrument must only be operated with a tube attached; with no tube there is a risk of injury from the rotating locating pin.

- Hearing protection must be worn when working with **BMT-20** or **BMT-50** tubes.

**DANGER**

Wear your personal protective equipment in accordance with the hazard category of the medium to be processed.

Otherwise there is a risk of:

- splashing liquids
- projectile parts
- body parts, hair, clothing and jewellery getting caught.

- Set up the instrument in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- The feet of the instrument must be clean and undamaged.
- The instrument is not suitable for manual operation.
- The instrument may heat up when in use.
- Check the instrument and accessories beforehand for damage

- each time when you use them. Do not use damaged components.
- Ensure that the cover on the tube is screwed on tightly.
- Ensure that the tube is firmly attached to the bayonet lock connector on the drive unit prior to operating the instrument.
- The tube must only be attached and removed while the motor is stationary.

**WARNING**

The **IKA**® tubes must always be closed when the instrument is in operation. Switch off the instrument immediately if any material leaks from the tube. Clean the instrument.

- Always open the tube carefully after use as the media in the tube may heat up due to transfer of energy during operation, leading to pressurization of the container: risk of material spraying, protective equipment must be worn.
- The temperature of the material must not exceed 40 °C.
- Only use tubes approved by **IKA**®.
- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
- Beware of the risk of flammable materials.
- **Do not** operate the instrument in explosive atmospheres, with hazardous substances or under water.
- The instrument doesn't start up again automatically following a cut in the power supply.
- Safe operation is only guaranteed with the accessories described in the "**Accessories**" chapter.

For protection of the instrument

- The voltage stated on the name plate must correspond to the mains voltage.
- The instrument must only be operated with the original plug-in power supply unit.
- Protect the instrument and accessories from bumps and impacts.
- The instrument may only be opened by experts.

## 4 Unpacking

### • Unpack

- Please unpack the device carefully
- In the case of any damage a fact report must be set immediately (for post, rail or forwarder)

### • Delivery scope


- **Tube Drive control**
- Power supply unit with exchangeable adapters
- Removal hook
- Operating instructions
- A warranty card
- A decontamination clearance certificate.

Power supply unit with exchangeable adapters.

Power supply

Exchangeable adapter  
Europe, Switzerland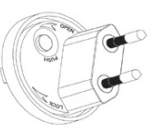

Exchangeable adapter  
USA, China

Exchangeable adapter  
England

Exchangeable adapter  
Australia

Optionally in the storage case:

- 1 x ST 20
- 1 x DT 20
- 1 x BMT 20 S
- 1 x BMT 20 G

## 5 Correct use

### • Use

	Application	Volume	Mode
▶ with Tube ST-2	Mixing		Seconds and minutes mode
▶ with Tube ST-20...	Stirring	1 - 2 ml 2 - 15 ml	
▶ with Tube ST-50...		15 - 50 ml	Seconds and minutes mode
▶ with Tube DT-20...	Dispersing		
▶ with Tube DT-50...	Stirring	5 - 15 ml 15 - 50 ml	minutes mode
▶ with Tube BMT-20...	Crushing		Seconds and minutes mode
▶ with Tube BMT-50...	Mixing we/dry	2 - 15 ml 15 - 50 ml	

### • Range of use (indoor use only)

- Laboratories
- Schools
- Pharmacies
- Universities

This device is suitable for use in all areas except:

- Residential areas
- Areas that are connected directly to a low-voltage supply network which also supplies residential areas.

The safety of the user cannot be guaranteed:

- If the instrument is operated with accessories that are not supplied or recommended by the manufacturer
- If the instrument is operated improperly contrary to the manufacturer's specifications
- If the instrument is operated improperly contrary to the manufacturer's specifications.


## 6 Useful facts

Changes in viscosity and volume caused by processing of dispersion media may cause small fluctuations in the speed of the device.


## 7 Commissioning

Pay attention to the ambient conditions listed in the “**Technical Data**”.

Position the rotating knob (B) at the left stop before starting up.

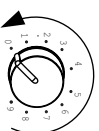
If these conditions are met, the unit is ready for service when the mains plug has been plugged in.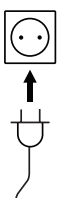

Fit the tube carefully and ensure it is properly attached.

If above procedures are not followed, safe operation can not be guaranteed and/or the equipment may be damaged.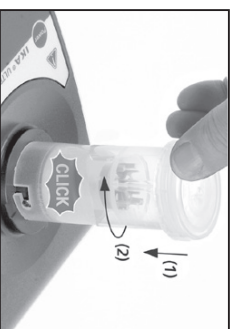


## Switching on

Press “Power” button (A).

Previous mode and time settings will be displayed e.g.:

(Factory setting: -2)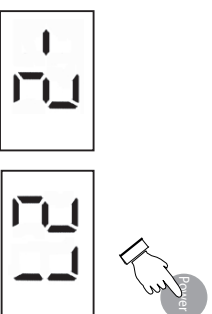


## Setting Timer (Changing of mod)

Switch of the device.

### Minutes-mode

Press and hold “Start/STOP” button (C), and also press the “Power” button (A).

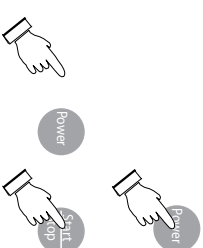Minutes-mode display:

The time that was previously set will be stored, e.g.:

*Note: max. 29 min., speed range upto 4000 min<sup>-1</sup>.*

### Seconds-mode

Press and hold “Start/STOP” button (C), and also press the “Power” button (A).

Seconds-mode display:

The time that was previously set will be stored, e.g.:


*Note: max. 59 sec, speed range upto 6000 min<sup>-1</sup>.*

## Setting the time

The time that was previously set will be stored, e.g.:

(Factory setting: 00)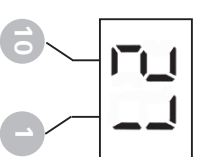


Set the tens digit using button (E).      Set the ones digit using button (D).

### Attaching the tube

Fit the tube carefully and ensure it is properly attached.

### Setting the speed

Set the desired speed in the range of 300 - 6000 min<sup>-1</sup> using the infinitely adjustable rotary knob (B).

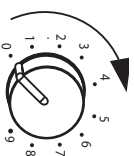### Starten

Press "Start/Stop" button (C), while the preset countdown is running, timer display will blink.

An audible signal will be given when the countdown is complete.

### Cancelling the countdown

Press button "Start/Stop" (C).

### Ausschalten

Press button "Power" (A).

## 8 Error message

It is possible for the media to become lodged between the cutters of the DT-tubes. The unit will automatically change the direction of rotation up to three times in an attempt to free the cutters. The unit will switch off automatically if this is not possible.

If the device becomes blocked on three occasions within a certain period of time the following error message will occur:

- ⇒ Display:
- ⇒ An audible signal will be given.

**Corrective action:** Switch off the device, check contents, switch on the device.


## 9 Accessories

- **ST-2**  
Tube with stirring device 2 ml
- **ST-20**  
Tube with stirring device 20 ml, with pierceable membrane
- **ST-20-M-gamma**  
Tube with stirring device 20 ml, γ-sterilized, with pierceable membrane
- **ST-50**  
Tube with stirring device 50 ml
- **ST-50-M**  
Tube with stirring device 50 ml, with pierceable membrane
- **ST-50-M-gamma**  
Tube with stirring device 50 ml, γ-sterilized, with pierceable membrane
- **DT-20**  
Tube with dispersing device 20 ml with pierceable membrane
- **DT-20 eco**  
Tube with dispersing device 20 ml, with PSU rotor/stator
- **DT-20-M-gamma**  
Tube with dispersing device 20 ml, γ-sterilized, with pierceable membrane
- **DT-50**  
Tube with dispersing device 50 ml
- **DT-50 eco**  
Tube with dispersing device 50 ml, with PSU rotor/stator
- **DT-50-M**  
Tube with dispersing device 50 ml, with pierceable membrane
- **DT-50-M-gamma**  
Tube with dispersing device 50 ml, γ-sterilized, with pierceable membrane

- **BMT-20-S** Tube for grinding with stainless steel balls 20 ml, with pierceable membrane
- **BMT-20-S-M-gamma** Tube for grinding with stainless steel balls, 20 ml, γ-sterilized, with pierceable membrane
- **BMT-50-S** Tube for grinding with stainless steel balls 50 ml
- **BMT-50-S-M** Tube for grinding with stainless steel balls 50 ml, with pierceable membrane
- **BMT-50-S-M-gamma** Tube for grinding with stainless steel balls 50 ml, γ-sterilized, with pierceable membrane
- **BMT-20-G** Tube for grinding with glass balls 20 ml, with pierceable membrane
- **BMT-50-G** Tube for grinding with glass balls 50 ml
- **BMT-50-G-M** Tube for grinding with glass balls 50 ml, with pierceable membrane
- **ST-2-adapter** Adapter UTTD - ST-2
- **TC-20-M** Tube cover 20 ml with pierceable membrane
- **TC-50** Tube cover 50 ml
- **TC-50-M** Tube cover 50 ml with pierceable membrane
- Glass balls
- Stainless steel balls

## 10 Maintenance

The drive unit is maintenance-free but not immune from wear. The carbon brushes of the motor wear down over the time.

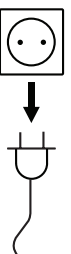For cleaning disconnect the main plug.

Only use cleansing agents which have been recommended by **IKA®**.

### Dirt

Dyes

Construction materials

Cosmetics

Foodstuffs

Fuels

### Cleaning agent

isopropyl alcohol

water containing tenside/isopropyl alcohol

water containing tenside/isopropyl alcohol

water containing tenside

water containing tenside

For materials which are not listed, please request information from **IKA®** application support.

Wear protective gloves during cleaning the instruments.

Electrical instruments may not be placed in the cleansing agent for the purpose of cleaning.

Do not allow moisture to get into the instrument when cleaning.

Before using another than the recommended method for cleaning or decontamination, the user must ascertain with **IKA®** that this method does not destroy the instrument.

### Spare parts order

When ordering spare parts, please give:

- machine type
- manufacturing number, see type plate
- item and designation of the spare part, see [www.ika.com](http://www.ika.com), spare parts diagram and spare parts list.

## Repair

**In case of repair the device has to be cleaned and free from any materials which may constitute a health hazard.**

For repair, use the form supplied or printed copy on the **IKA®** website: **www.ika.com**.

If you require servicing, return the appliance in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

## 11 Warranty

In accordance with **IKA®** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

## 12 Technical data

### Power supply unit

Input	<b>VAC</b> 100 - 240
	<b>A</b> 0,6
	<b>Hz</b> 50-60
Output	<b>VDC</b> 24
	24W,LPS (limited power source)
Protection class	2 (double insulated) 

### Tube Drive

Operating voltage	<b>VDC</b> 24
	<b>mA</b> 1000
Power consumption, normal operation	<b>W</b> 20
Power consumption, standby operation	<b>W</b> 2
Power output	<b>W</b> 17
Drive	EC flat motor
Speed range	<b>rpm</b> 400 - 6000 infinitely adjustable, adjustment magnitude 10 rpm up to 8000
Turbo speed	up to 8000
Mode A	up to 6000
Mode B	up to 6000
Mode P1-P9	up to 6000
Speed setting	rotating/pressing knob
Speed display	digital
Timer	10 seconds-30 minutes
Reverse timer	10 - 60 Seconds
Display-timer	digital
Perm. ambient temperature	<b>°C</b> +5 to +40
Perm. rel. humidity	<b>%</b> 80
Protection to DIN EN 60529	IP 20
Contamination level	2
Overvoltage category	II
Operation at a terrestrial altitude level	<b>m</b> max. 2000
Dimensions (W x D x H) without tube	<b>mm</b> 122 x 178 x 55
Weight	<b>kg</b> 1,0

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**IKA® - Werke GmbH & Co.KG**