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- Read the operating instructions in full before starting up and follow the safety instructions.**
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- CAUTION!** For research purposes only. Not suitable for therapeutic or diagnostic use in human or veterinary medicine (in accordance with the IVD Directive).
- Ensure that the cover is screwed on tightly.
- The tube must only be attached and removed whilst the motor is stationary.
- Ensure that the tube is firmly attached to the bypass lock connector on the drive unit prior to operating the device.
- Ensure that the tube is able to withstand the chemical properties of the selected medium by filling the tube and stirring for a suitable period time.
- Always open the tube carefully after use as the media in the tube may heat up due to transfer of energy during operation, leading to pressurization of the container, risk of material spraying, protective equipment must be worn.
- Abrasion of the dispersion equipment or the rotating accessories can get into the medium you are working on.
- While working with the dispenser, the user must wear his personal protective equipment, in particular safety goggles.
- The temperature of the material must not exceed 40 °C.
- Note the operating instructions of the dispenser unit.

Product information
<p>The operation of the ST1-2 on the dispenser is only possible with the ST1-2 adapter. A ST1-2 adapter is included. ST1/ST1-20/50... comprise mixing vessel and stirring tool, DT/20/50... comprise mixing vessel and dispersion tool.</p> <p>BMT/DT/5050... comprise mixing vessel, stirring tool and balls.</p> <p>ST1-20... M1-50... ST1-50-M... DT-20... DT-20-M... DT-50-M... DT-50-M and BMT-20... M1-50... BMT-50-M... BMT-50-M are not sterile and not free of DNA or RNA.</p> <p>ST1-20-gamma, ST1-20-M-gamma, DT-20-gamma, DT-20-M-gamma and BMT-20-gamma, BMT-20-M-gamma are sterilized.</p> <p>TC-2-M1, TC-20-M1 and TC-50-M are tube covers with a permeable membran for needles or pipette points.</p> <p>The tubes are ideal for applications, with which no cross contamination is permissible.</p> <p>The tubes ST1-20, DT1-20 and BMT1-20 are appropriate only for a single use. Multiple use can destroy the tube and the dispenser unit.</p>
Application instruction

Store the tubes in a dry place, in their packaging and at room temperature.

Note the use-by date of the tubes.

Protect the tubes against ultraviolet light.

Attention! Never run the DT-20/50... dry.

For ST1-2 the fill level of the medium must be between the 1 ml and 2 ml marks, for ST 20.../DT-20... and BMT-20... the fill level of the medium must be between the 2 l, respectively 5 ml and max. 15 ml marks, for ST 50.../DT-50... and BMT-5050... the fill level of the medium must be between the 15 ml and max. 50 ml marks.

Don't turn the tube after piercing the membran, then the leak-tightness is ensured no longer.

Always position the sample in the centre of the dispersing unit.

The solid bodies in samples must not be frozen or too hard.

Dispensing causes the medium to heat up.

Use the optimal dispersion, duration and rotating frequency must be determined by attempts. Usually a few seconds are sufficient.

Longer operation is not recommended.

Further application brings no improvements, increase however the sample temperature substantially.

Dispose of the tubes after use according to the regulations. Observe the laboratory regulations.

Materials and chemical resistance
<p>The containers are manufactured from polypropylene (PP), polyethylenketone (PEEK), Rotorsatzor are manufactured from polyethylenketone (PEEK) and polyethylenketone with teflon (PEEK/PTFE). At DT/20/50 eco are rotorsatzor of polyethersulfone (PESE) and polyethylenketone with teflon (PEEK/PTFE).</p> <p>The balls are manufactured from either stainless steel AISI 304 or soda-line glass.</p> <p>The used plastic materials conform with the FDA, food and Drug Administration Approval)... rules.</p> <p>The parts have a good stability against weak acids, chlorides, hypochlorides and many other chemicals.</p>
Technical data

Designation	For dispenser unit	Working range	Stator ø	Rotor ø	Circumferential speed at 6.000 rpm
ST1-2	Tube Drive	1-2 ml	-	-	-
ST1-20...	Tube Drive	2-15 ml	-	-	-
DT1-20...	Tube Drive	5-15 ml	26 mm	20 mm	6,5 m/s
BMT-20...	Tube Drive	2-15 ml	-	-	-
ST1-50...	Tube Drive	15-50 ml	-	-	-
DT1-50...	Tube Drive	15-50 ml	26 mm	20 mm	6,5 m/s
BMT-50...	Tube Drive	15-50 ml	-	-	-

Warranty

In accordance with **KA[®]** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer, you may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

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