**IKA® Vortex Genius 3**

Betriebsanleitung Ursprungssprache	DE	03
Operating instructions	EN	07
Mode d'emploi	FR	11
Руководство пользователя	RU	15

Indicaciones de seguridad	ES	<b>19</b>	Biztonsági utasítások	HU	<b>29</b>
Veiligheidsinstructies	NL	<b>20</b>	Varnostna opozorila	SL	<b>30</b>
Norme di sicurezza	IT	<b>21</b>	Bezpečnostné pokyny	SK	<b>31</b>
Säkerhetsanvisningar	SV	<b>22</b>	Ohutusjuhised	ET	<b>32</b>
Sikkerhedshenvisninger	DA	<b>23</b>	Drošības norādes	LV	<b>33</b>
Sikkerhetsanvisninger	NO	<b>24</b>	Nurodymai dėl saugumo	LT	<b>34</b>
Turvallisuusohjeet	FI	<b>25</b>	Инструкции за безопасност	BG	<b>35</b>
Instruções de serviço	PT	<b>26</b>	Indicații de siguranță	RO	<b>36</b>
Wskazówki bezpieczeñstwa	PL	<b>27</b>	Υποδειξείς ασφάλειας	EL	<b>37</b>
Bezpečnostní pokyny	CS	<b>28</b>			


Reg.-No. 4343-01

## Bedienelemente / Control elements / Éléments de réglage / Коммутационные элементы

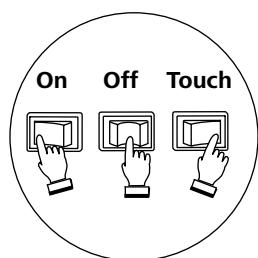Fig. 1

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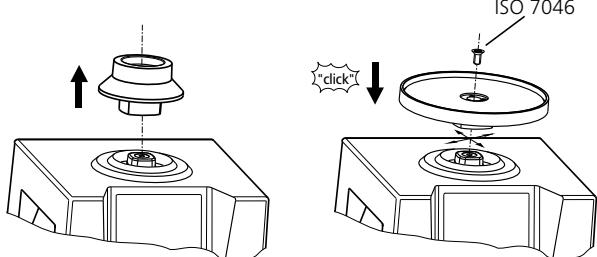
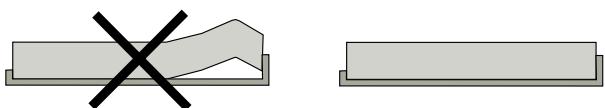
Fig. 4

Fig. 3

- Reduce the speed if
- the medium splashes out of the vessel because the speed is too high
- the appliance is not running smoothly
- the appliance begins to move around because of dynamic forces.

#### Accessories

- Safe operation is only guaranteed with the accessories described in the "Accessories" chapter.

- Always disconnect the plug before attaching accessories.
- Accessories must be securely attached to the device and cannot come off by themselves. The centre of gravity of the assembly must lie within the surface on which it is set up.

#### For protection of the equipment

- The appliance may only be opened by experts.
- Protect the appliance and accessories from bumps and impacts.

## Unpacking

### • Unpacking

- Please unpack the device carefully
- In the case of any damage a report must be sent immediately (post, rail or forwarder).

### • Scope of delivery

- Shaker **Vortex Genius 3**
- Standard attachment VG 3.1
- Operating instruction

## Correct use

### • Use

- For mixing liquids
- Touch mode for single test tubes
- Continuous mode for single or several vessels

### • Range of use (only indoors)

- Laboratories
- Schools
- Pharmacies

This device is suitable for use in all areas except:

- Residential areas
- Areas that are connected directly to a low-voltage supply network that also supplies residential areas.

The safety of the user cannot be guaranteed if: the appliance is operated with accessories that are not supplied or recommended by the manufacturer, the appliance is operated improperly according to the manufacturer's specifications or the appliance or printed circuit board are modified by third parties.

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**IKA® - Werke GmbH & Co.KG**