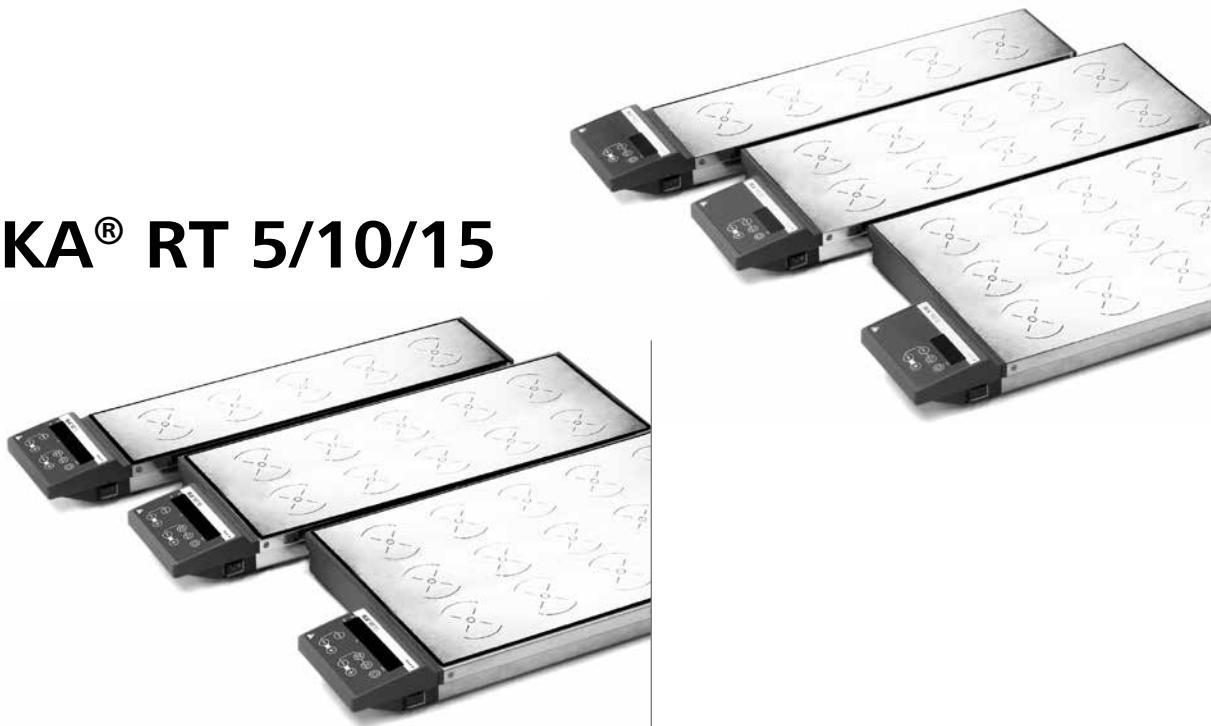
IKA® RO 5/10/15**IKA® RT 5/10/15**

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	Page
Control panel and display	02
Contents	09
Declaration of conformity	09
Warranty	09
Explication of warning symbols	09
Safety instructions	10
Unpacking	11
Correct use	11
Commissioning	11
Safe temperature limit	11
Setting the safe temperature limit	11
Stirring function	11
Eco mode	12
Operating modes	12
Direction of rotation reversal	12
Function heating	12
Maintenance and cleaning	12
Error codes (RT 5/10/15)	13
Accessories	13
Technical data	14

Declaration of conformity

We declare under our sole responsibility that this product is in compliance with the regulations 2006/95/EC and 2004/108/EC and conforms to the standards or standardized documents
DIN EN IEC 61010-1, -2-010, -2-051 and DIN EN IEC 61326-1.

Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

Explication of warning symbols

General hazard.

This symbol identifies information that is of absolute importance to ensure health and safety. Failure to observe this information may be detrimental to health or may result in injuries.

This symbol indicates information which is important for ensuring that the appliance functions without any technical problems. Failure to observe this information could damage the appliance.

This symbol indicates information which is important for proper use of the appliance and / or ensuring that the appliance functions correctly. Failure to observe this information can lead to inaccurate results.

ATTENTION - Risk of damage due to magnetism.

DANGER - Reference to the endangerment by a hot surface

Safety instructions

For your protection

- **Read the operating instructions in full before starting up and follow the safety instructions.**
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Socket must be earthed (protective ground contact).

• **Attention - Magnetism!**

Effects of the magnetic field have to be taken into account (e.g. data storage media, cardiac pacemakers ...).

RT

• **Risk of burns!**

Exercise caution when touching the housing parts and the heating plate.

The heating plate can reach temperatures in excess of 125 °C. Pay attention to the residual heat after switching off.

RT

- Ensure that the mains power supply cable does not touch the heating base plate.

- Wear your personal protective equipment in accordance with the hazard category of the media to be processed. Otherwise there is a risk from:
 - splashing and evaporation of liquids
 - ejection of parts
 - release of toxic or combustable gases.
- Set up the appliance in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- The feet of the appliance must be clean and undamaged.
- Check the appliance and accessories beforehand for damage each time you use them. Do not use damaged components.
- Gradually increase the speed.
- Reduce the speed if
 - the medium splashes out of the vessel because the speed is too high
 - the appliance is not running smoothly
 - the container moves on the base plate.

RT

- **Caution!** Only process and heat up any media that has a flash point higher than 180°C.

The safe temperature limit must always be set to at least 25 °C lower than the fire point of the media used.

- Beware of hazards due to:

- flammable materials
- combustible media with a low boiling temperature
- glass breakage
- incorrect container size
- overfilling of media
- unsafe condition of container.
- The appliance may heat up when in use.
- The base plate can heat up due to the action of the drive magnets at high motor speeds, even if the heater is not operational.

- Process pathogenic materials only in closed vessels under a suitable extractor hood. Please contact **IKA** if you have any questions.

DANGER

DANGER

plies to any extra energy produced in other ways, e.g. through light irradiation.

- Please observe the operating instructions for any accessories used.
- Safe operation is only guaranteed with the accessories described in the „Accessories“ chapter.
- Accessories must be securely attached to the device and cannot come off by themselves. The centre of gravity of the assembly must lie within the surface on which it is set up.
- Always disconnect the plug before fitting accessories.
- The appliance can only be disconnected from the mains supply by pulling out the mains plug or the connector plug.
- The socket for the mains cord must be easily accessible.
- It may be possible for wear debris from rotating accessory parts to reach the material being processed.
- When using PTFE-coated magnetic bars, the following has to be noted:

Chemical reactions of PTFE occur in contact with molten or solute alkali metals and alkaline earth metals, as well as with fine powders of metals in groups 2 and 3 of the periodic system at temperatures above 300 °C - 400 °C. Only elementary fluorine, chlorotrifluoride and alkalinmetals attack it; halogenated hydrocarbons have a reversible swelling effect.

(Source: Römpps Chemie-Lexikon and „Ullmann“ Volume 19)

For protection of the equipment

- The appliance may only be opened by experts.
- The voltage stated on the type plate must correspond to the mains voltage.
- Do not cover the device, even partially e.g. with metallic plates or film. This results in overheating.
- Protect the appliance and accessories from bumps and impacts.
- Ensure that the base plate is kept clean.
- Observe the minimum distances between the devices, between device and wall and minimum distances (min. 800 mm) above the assembly

see Fig. 2.

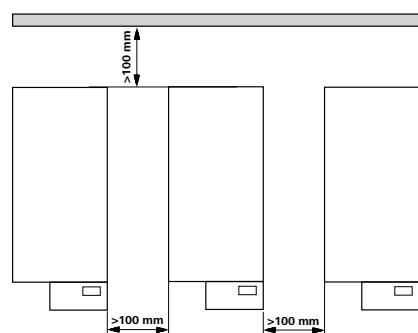
Fig. 2

Unpacking

• Unpacking

- Please unpack the device carefully
- In the case of any damage a detailed report must be set immediately (post, rail or forwarder)

• Delivery scope

RO

- Magnetic stirrer RO5/10 or 15
- power supply unit
- operating instruction

RT

- Heating magnetic-stirrer RT5/10 or 15
- mains cable
- operating instruction

Correct use

• Use

- For mixing and / or heating liquids

• Range of use (inside only)

- | | |
|----------------|----------------|
| - Laboratories | - Schools |
| - Pharmacies | - Universities |

The device is suitable for use in all areas except:

- Residential areas

- Areas that are connected directly to a low-voltage supply network that also supplies residential areas.

The safety of the user cannot be guaranteed if the appliance is operated with accessories that are not supplied or recommended by the manufacturer or if the appliance is operated improperly contrary to the manufacturer's specifications or if the appliance or the printed circuit board are modified parties.

Commissioning

RO/RT

Observe the ambient conditions (temperature, humidity, etc.) listed under Technical Data.

The device is switched on and off via the switch (A) located on the right-hand side of the device. On the front panel a small arrow labelled Power above the display (B) indicates the location of the switch.

After the device is switched on a display self-test is performed. The display cycles through the following information:

- All segments light up
- The software version and the selected mode are displayed
- The selected "Eco Mode" is displayed
(E when ECO mode is active)
- If the drive has not been started the display reads „OFF". If the drive has been started the selected speed is displayed.

Safe temperature limit

RT

The max. attainable heating plate temperature is restricted to 130°C by a permanently set safety temperature limit. Once this limit has been attained, the device switches off the heating.

WARNING
The safe temperature limit must always be set at least 25 °C lower than the flash point of the media to be processed!

Setting the safe temperature limit

RT

The safety temperature is set as follows:

- Switch the device on using the On/Off switch on the right side of the device.
- As soon as the display shows SAFE, press and hold down the Temp membrane key and use

the +/- membrane keys (**H** or **I**) to set the desired safety temperature.

The set safety temperature is displayed.

Once the Temp membrane key is released, the safety temperature is saved.

Stirring function

RO/RT

The stirring function is started and stopped by pressing the „Start/Stop" button (**C**).

Pressing the +/- buttons (**E or D**) increases or reduces the speed.

ECO mode

RO/RT

The heat generated in the device can be reduced by lowering the power of the input coils.

Pressing the „ECO Mode“ (**F**) key reduces the current power setting. ECO mode is indicated by an “E” preceding the speed. When the power is set to ECO mode, the speed is limited to 600 rpm.

Operating modes

RO/RT

To toggle between modes, press the Start/Stop key (C) when switching on.

Operating Mode “A”

All the parameter settings are retained when the device is switched off or disconnected from the mains.

The “Start/Stop” (**C**) key starts the device with the set parameters.

Operating Mode “B”

All parameter settings, including “Start/Stop”, are saved when the device is switched off.

Factory setting: mode A

Direction of rotation reversal

RO/RT

The (**G**) key activates the “Reverse Rotation” function. When the „Reverse Rotation“ function is switched on, a decimal point (**K**) appears in the LCD display.

The running time (30, 60, 90 or 120sec) is set by pressing the (**G**) key followed by the +/- keys.

After the preset running time, the drive is switched off and restarted at minimum speed in the opposite direction after a pause specific to the speed.

In this way the magnetic rod is captured and any liquid still rotating is gradually slowed down. Then the drive accelerates to the target speed.

Note: If the “Reverse Rotation” key (**G**) is pressed again, the drive will keep rotating in the same direction.

When the device is started up again, the drive will rotate in the direction last set.

Function Heating

RT

The heating function is started and stopped by pressing the „Temp“ button (**J**).

Pressing the +/- buttons (**H** or **I**) increases or reduces the temperature.

Maintenance and cleaning

The equipment is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.

Cleaning

For cleaning disconnect the main plug..

Use only cleansing agents which have been approved by **IKA** to clean devices.

Use to remove:

Dyes	Isopropyl alcohol
Construction materials	water containing ten side / isopropyl alcohol
Cosmetics	water containing ten side / isopropyl alcohol
Foodstuffs	water containing tenside
Fuels	water containing tenside
- Do not allow moisture to get into the appliance when cleaning.	
- Wear protective gloves when cleaning the devices.	
- Please consult with IKA before using any cleaning or decontamination methods, other than those recommended here.	

Ordering spare parts

When ordering spare parts, please give:

- Device type
- Manufacturing number, see type plate
- Software version
- Item number and designation of the spare part,
see www.ika.com.

Repair

The device must be clean and free from any materials which may constitute a health hazard when sent for repair.

Use the „certificate of compliance“ included or download the form from the IKA® website www.ika.com.

Please return the appliance in its original packaging. Storage packaging is not sufficient for returns. Please also use suitable packaging for transportation.

Error Codes (RT 5/10/15)

Error-code	Cause	Effect	Solution
Er 3	Temperature inside device too high	Heating off	- Switch off device and allow to cool down
Er 6	Break in safety circuit	Heating off	- Plug in contact plug (L) - Plug in contact thermometer or PT 1000 temperature sensor - Replace faulty connection cable, plug, or contact thermometer
Er21	Safety relay doesn't open during test		- Switch off device <i>- Warning! Only to be carried out by authorised service personnel: Check PCB</i>
Er22	Safety temperature is too low during test	Heating off	- Switch off device <i>- Warning! Only to be carried out by authorised service personnel: Check PCB and safety temperature sensor</i>
Er25	Heating and switching element monitoring	Heating off	- Check the external temperature controller (f.e. Er 5 ETS-D) - Switch off device <i>- Warning! Only to be carried out by authorised service personnel: Carry out an internal test on the device to check the plug-in connector for the heating element, the safety temperature sensor or the PCB</i>
Er26	Difference between temperature of safety sensor and temperature of control sensor control temperature > (safety temperature + 40 K)	Heating off	- Switch off device <i>- Warning! Only to be carried out by authorised service personnel: Carry out an internal test on the device to check the plug-in connector for the temperature sensor</i>
Er31	Triac error detection doesn't work	Heating off	- Switch off device <i>- Warning! Only to be carried out by authorised service personnel: Check PCB</i>
Er46	Difference between temperature of safety sensor and temperature of control sensor control temperature > (control temperature + 40 K)	Heating off	- Switch off device <i>- Warning! Only to be carried out by authorised service personnel: Carry out an internal test on the device to check the plug-in connector for the temperature sensor</i>

If the actions described fail to resolve the fault or another error code is displayed then take one of the following steps:

- Please contact the service department;
- Send the device for repair, including a short description of the fault.

Accessories

- Stirring bar: ø 8mm; lenght 30 mm
- RSE Stirring bar remover

Technical data

RO

Power supply unit

Input	V	100 - 240	
Output	A	1,67	
	Vdc	24; 40W.LPS; (limited power source)	
Protection class	Hz	50/60	
Operating voltage		2 (double insulated) 	
	Vdc	24	
	RO 5	RO10	RO15
Power consumption (max.)	A	1,0	1,5
Power consumption, standby operation	W	24	36
Power output	W	2	
Speed ranges	W	17	
P100	rpm	0 - 1200 in 10rpm-steps adjustable	
ECO-Mode	rpm	0 - 600 in 10rpm-steps adjustable	
	RO 5	RO10	RO15
Set-up surface	mm	120 x 450	270 x 450
Dimension:	mm	120 x 570 x 60	280 x 570 x 60
Weight:	kg	3,0	7,0

RT

Unit

Operating voltage	Vac	230 ± 10%	
Frequency	Vac	115 ± 10%	
Speed ranges	Hz	50 / 60	
P100	rpm	0 - 1000 in 10rpm-steps adjustable	
ECO-Mode	rpm	0 - 600 in 10rpm-steps adjustable	
	RT 5	RT 10	RT 15
Power consumption (max.)	W	395	600
Set-up surface	mm	180 x 495	270 x 495
Dimension:	mm	190 x 610 x 60	280 x 610 x 60
Weight:	kg	4,0	9,4
Adjustment and display resolution	K	1	
Surface temperature	°C	120*	
Safety temperature limit (adjustable)	°C	50 - 150	
	RO/RT 5	RO/RT 10	RO/RT 15
Fuse:	230 V A	T 2 (5x20)	T 6,3 (5x20)
	100 V / 115 V A	T 4 (5x20)	T 10 (5x20)

RO/RT

Setting resolution	rpm	±5	
Deviation between the stirrer points	%	0	
Permissible duration of operation	%	100	
Permissible ambient temperature	°C	+5 to +40	
Permissible relative humidity	%	80	
EN 60529 protection class		IP 40	
Operation at a terrestrial altitude	m	max. 2000	
Stirred quantity max. (water)		400ml per stirrer point in the 600ml beaker glass	

	RO/RT 5	RO/RT10	RO/RT15
in other vessels	ltr	2	4
Stirrer points		5	10
Distance from stirrer point to stirrer point	mm	90	90 x 90

*Caution! The specified maximum hotplate temperature refers to instruments with a full load.
Without a load, the (setup) plate temperature can reach 130°C (+10°C / -5°C).
Load to determine the specified values: 250 ml Erlenmeyer flask filled with 200 ml water.

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IKA®-Werke GmbH & Co.KG