



Plastic Tray Assembly Instructions

PLEASE READ CAREFULLY: Proper assembly will ensure tray integrity and prolong life span.


1. Place the flat sheet so that the **scored grooved lines are facing up**. (Note: The scored grooved lines indent into the flat sheet)

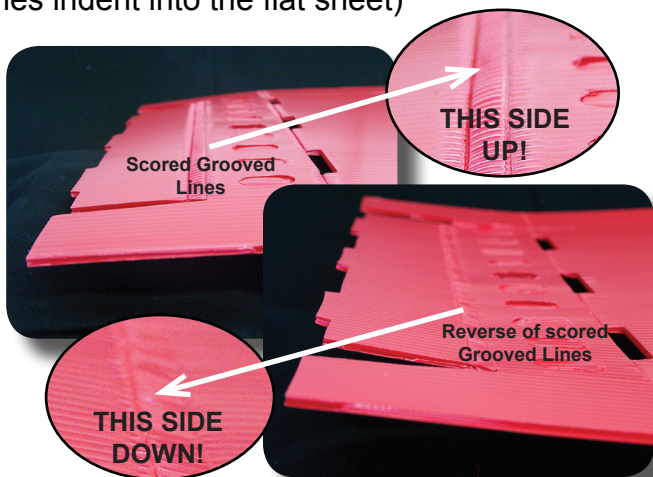


2. To facilitate folding the tray, pre-fold along all the scored grooved lines.

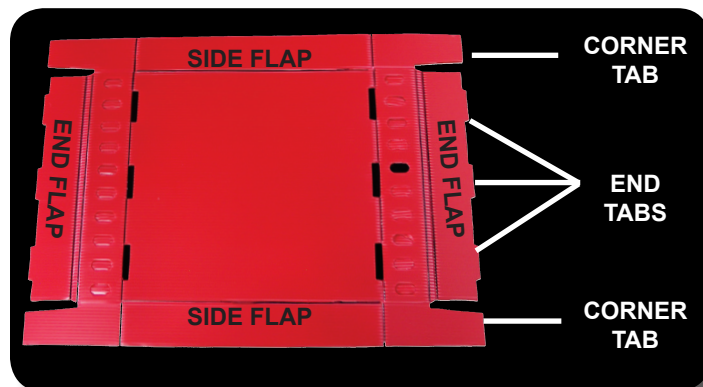


3. Lift the two side flaps up.

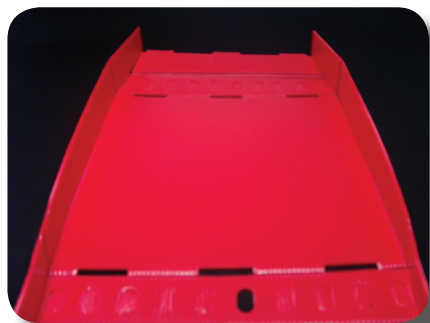

4. Fold the two corner tabs in so that they are perpendicular with the side flaps.

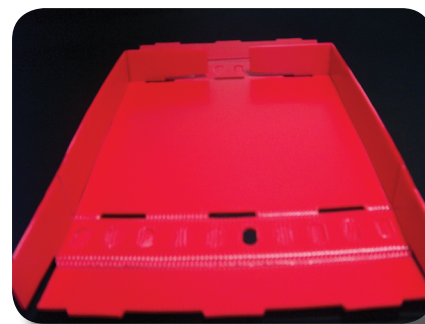

5. Lift the end flap up and over the corner tabs so that the end tabs fit into the holes on the bottom.

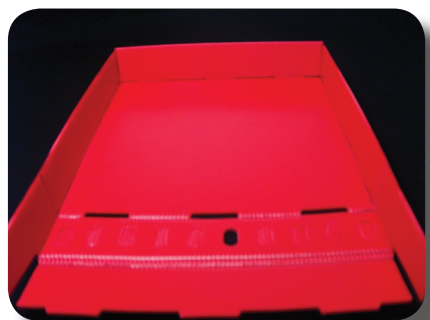

6. Repeat Steps 4 and 5 for the bottom wall.

